MTC695

User Manual

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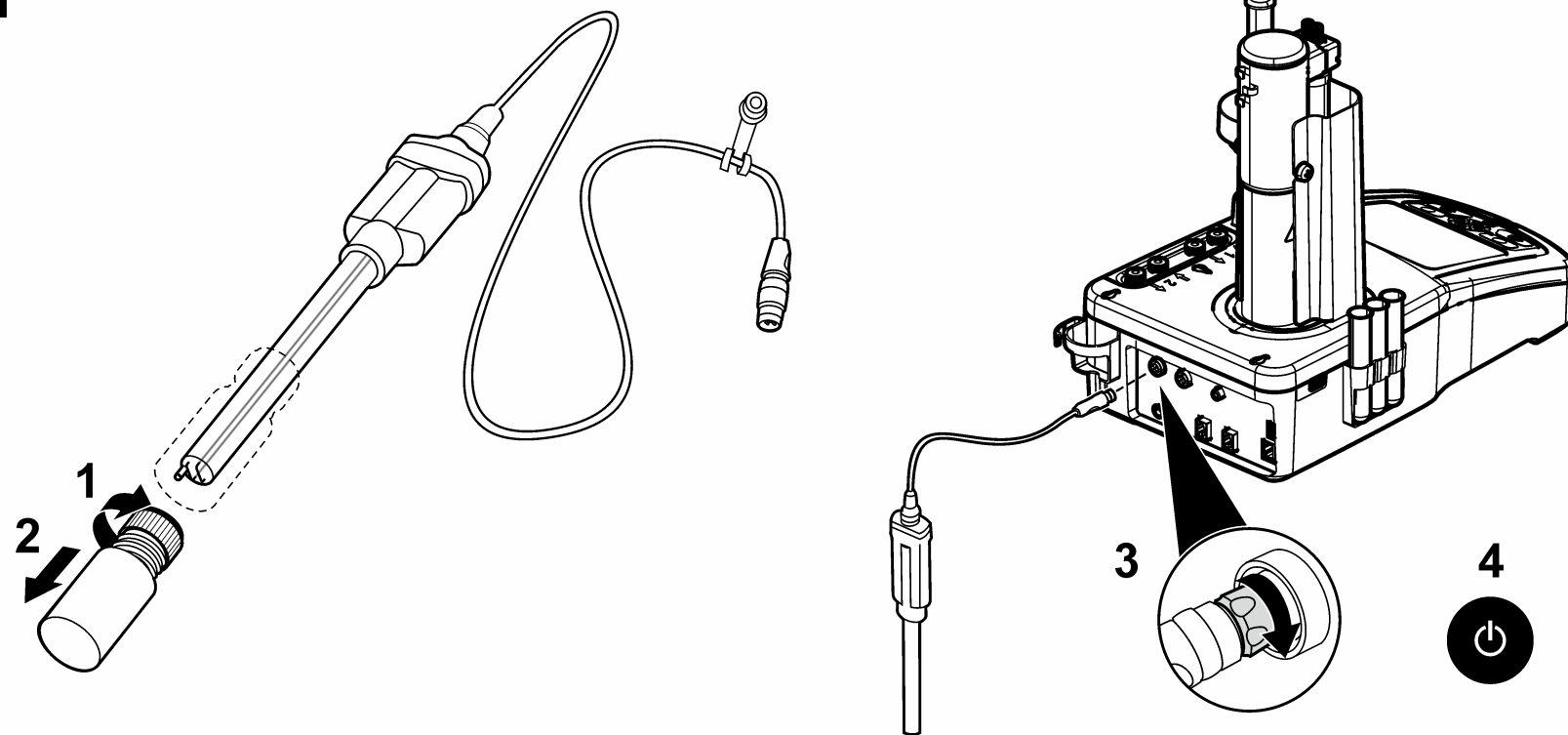
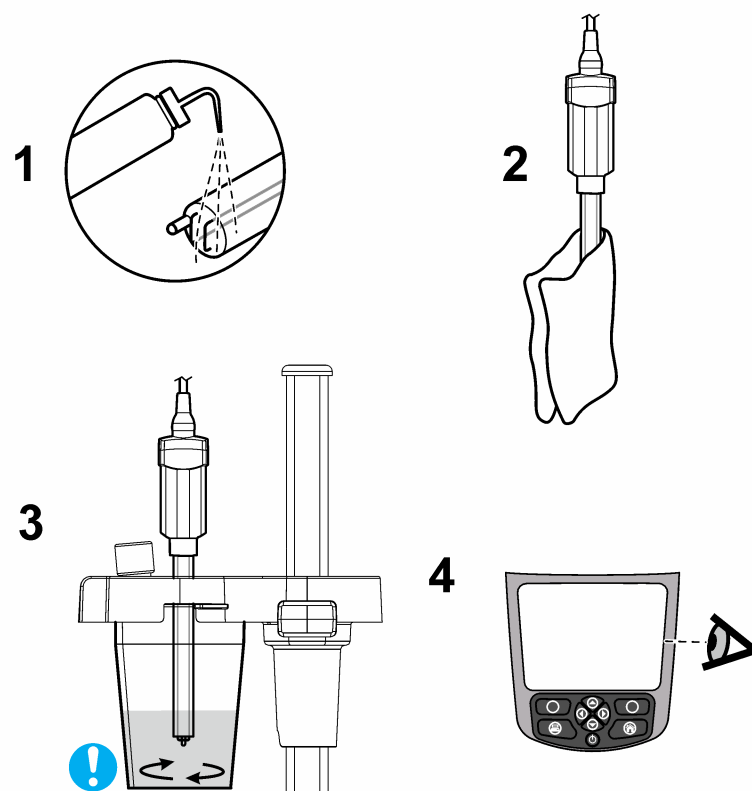
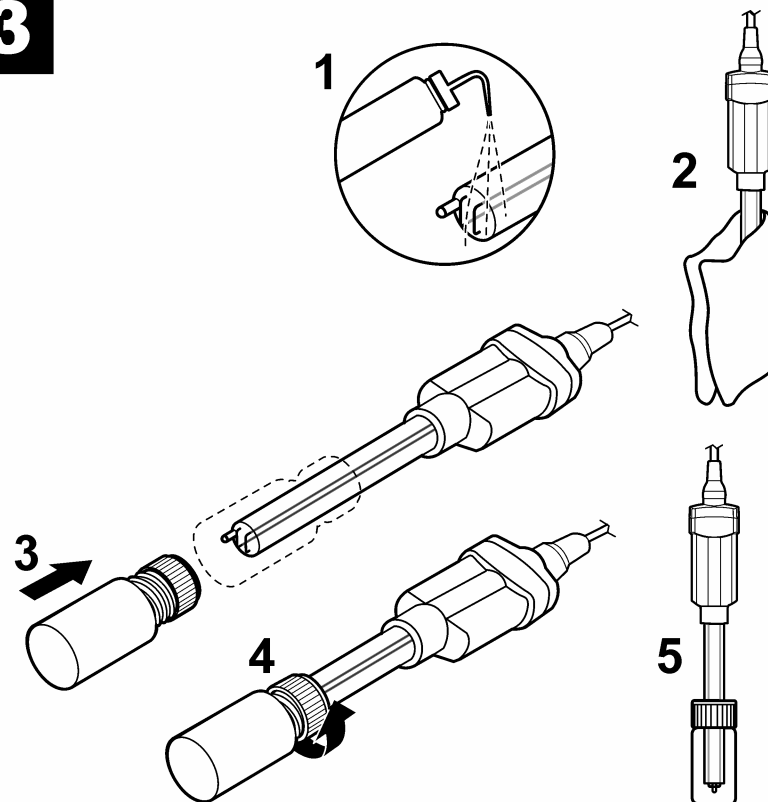
www.hach.com

1**2****3****EN****MTC695**

Additional information is available on the manufacturer's website.

Introduction

The probe is a double platinum electrode with a glass body and an integrated temperature sensor. The probe is used in Karl Fischer titrations and other potentiometric and amperometric titrations.

Warning

⚠ Chemical exposure hazard. Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.

Items to collect:

AT1000 or KF1000 series titrator, suitable solvent for probe cleaning, no-lint cloth

1 Connect the probe**2 Measurement****3 Storage**

For storage, keep the electrode dry in the soaker bottle at ambient temperature.

Refer to the instrument documentation for more information.

Technical information

Range	±1200 mV (±200 µA)
Temperature	-10 to 80 °C (14 to 176 °F)
Filling solution	-
Storage solution	None (dry)

Maintenance

Clean the probe with the applicable solvent solution. Refer to the applicable titration procedure (e.g., ethanol-based solvent for Karl Fischer titrations).

If necessary, use finishing abrasive bands to clean the platinum wires. Do not bend the platinum wires.

Warranty

1 year for manufacturing faults. This warranty does not cover improper use or wear.

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