ISE NH₄⁺ 9663C

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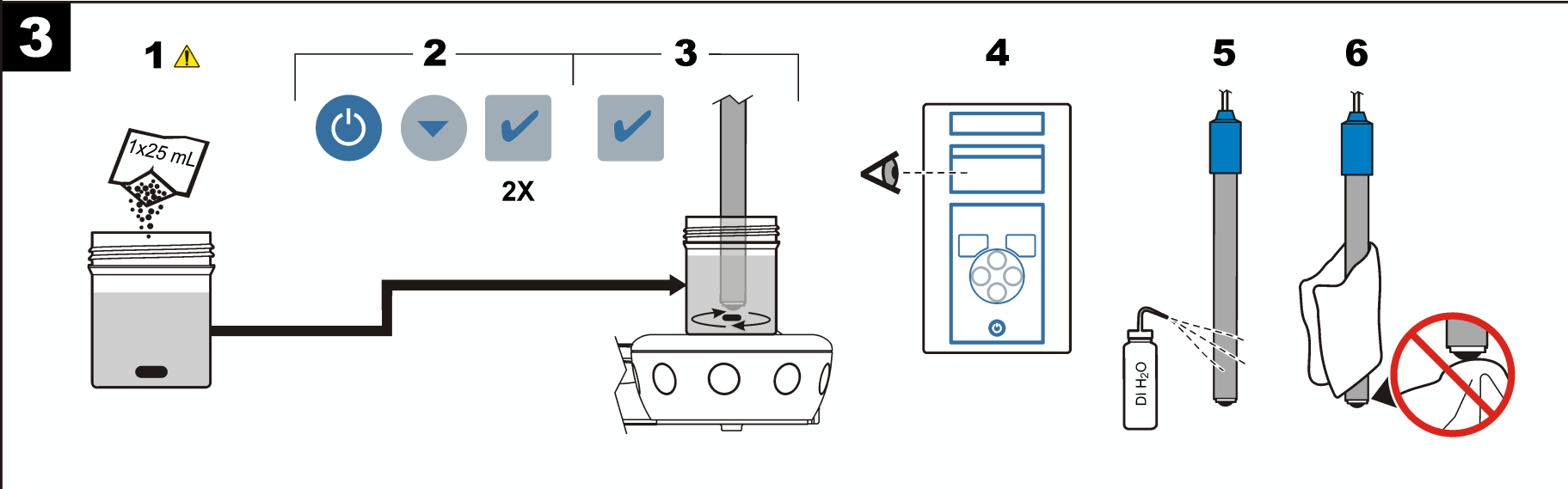
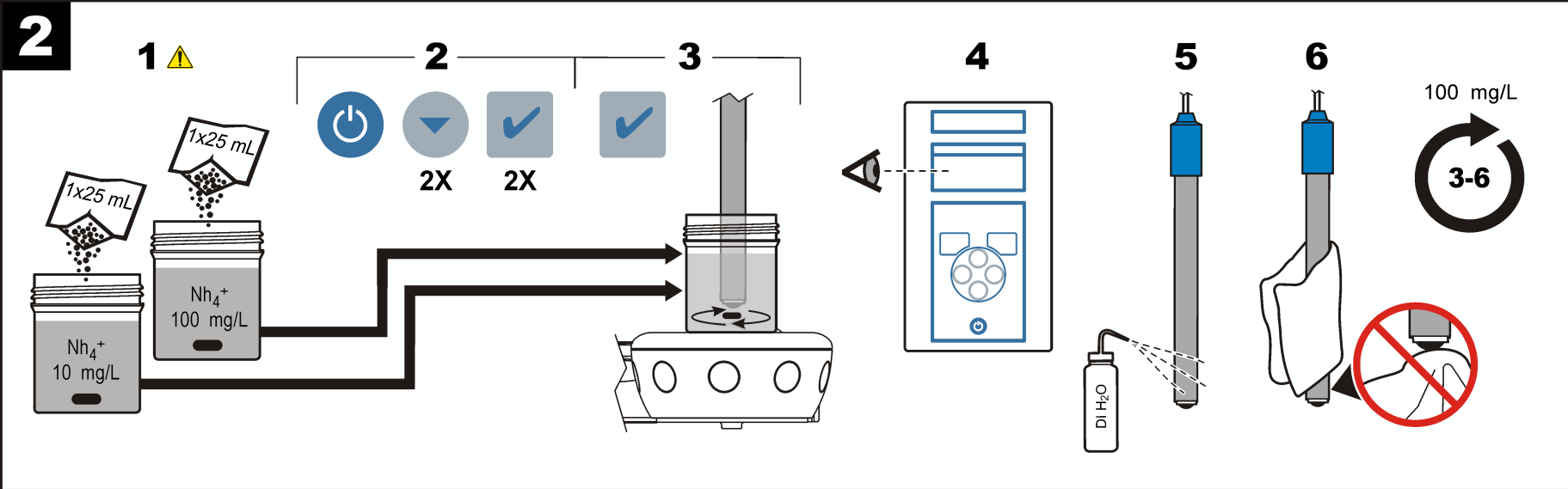
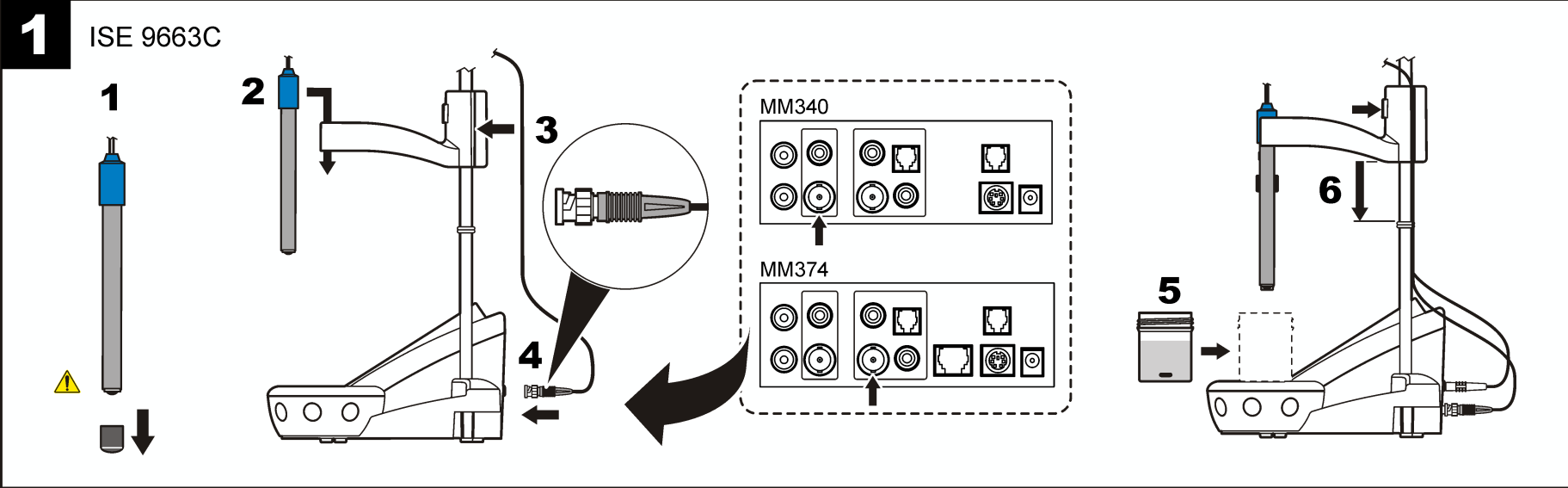
ISE NH₄⁺ 9663C

Refer to the meter manual for more information.

Caution

⚠ *Personal injury hazard. Glass components can break. Handle with care to prevent cuts.*

⚠ *Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.*

Introduction

This probe is a combination ammonium selective electrode with a non refillable double junction reference. The probe measures ammonium concentration in aqueous solutions.

Technical information

Range	0.018 mg/L (10 ⁻⁶ M) to 9,000 mg/L (0.5 M) as ammonium
pH range	0 to 8.5
Linear range	0.9 mg/L to 9,000 mg/L as Ammonium
Slope	58 mV/decade, 90 to 110% at 25 °C (77 °F)
Ionic strength adjuster (ISA)	Hach ISA 2980699
Operating temperature	5 to 50 °C (32 to 122 °F)
Interferences	K ⁺ (0.11), Na ⁺ (0.0016), Ca ²⁺ (0.00006), Li ⁺ (0.00003)

Preparation

1 Connect the probe.

2 Calibration

Add 1 ISA pillow per 25 mL of standard.

3 Measurement

Add 1 ISA pillow per 25 mL of sample.

Maintenance and storage

Cleaning: rinse with distilled water between each standard and each sample.

Storage: clean and dry the probe and install the sensor protection cap for storage.

Troubleshooting

The ionic strength and temperature of samples and standards must be the same. Make sure that the stir speed is the same for all standards and samples.

Damage or deposits on the membrane change the response of the electrode. Remove deposits from the membrane and replace the sensor if damaged.

Warranty

6 months for manufacturing faults only. The warranty does not cover normal wear or improper use.

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